

How to install Thin Client on IIS 6 / Windows Server 2003 32-bit

STEP 1:

Go to a command prompt.

Change to C:\inetpub\Adminscripts folder. (If Windows Server 2003 is installed on a drive other than "C:", use that drive letter instead.)

In the Adminscripts folder, type the following command:

```
cscript adsutil.vbs set w3svc/AppPools/Enable32bitAppOnWin64 1
```

Note: The value "1" for Enable32bitAppOnWin64 tells Windows to run the w3svc in 32-bit mode, while a value of "0" indicates 64-bits.

Type iisreset.

Reboot the server.

STEP 2:

ASP pages are not enabled in the default install of iis. Use add/remove programs to drill into the www publishing service config and select/install this functionality. Otherwise, none of the thin client will work.

STEP 3:

Run www service in IIS 5.0 isolation mode:

Open IIS manager

Open the Tree View

Right Click the "Web Sites" folder

Select Properties

Select the Services Tab

Click the "Run WWW Service in IIS 5.0 isolation mode" checkbox

Click Apply

Restart IIS if it does not automatically (reboot the server if you are not sure how to restart IIS)