

DV3111F – Report: System Configuration

Scope The System Configuration Report shows the configuration of every input on the system grouped by each Asset on the system.

Purpose The report can be used to check multiple Assets at a time to make sure that they are configured properly. See related documents for how to set up and edit Asset configuration.

Applicable User Groups All Users (Administrator, Group Administrator, General User)

Requirements User must be logged into the system and assigned to a group that has Assets assigned to the group as well.

The report includes the following fields:

1. Asset Name (Tag) (customizable asset name and unique asset tag)
2. Type (the type of appliance being monitored)
3. Location (Asset location)
4. Groups (list of the groups the Asset is assigned to)
5. Status (Active/Inactive)
6. Alarm Notification Schedule (name of the assigned alarm notification schedule)
7. Holiday Alarm Notification Schedule (name of the assigned holiday alarm notification schedule)
8. Input (information on the specific input including type of readings, input channel, device code, etc.)
9. Alarm Settings (including Min, Max, Alarm Delays and Latching settings for Pre-Alarm, Standard Alarm, and Emergency Alarm thresholds)

System Configuration						
Assets						
Asset Name (Tag)	Type	Location	Groups	Status		
AA04872 - Conf Rm (Ref) (AA04872)	Reach-In	Conf Room	global and Client Security	Active		
Alarm Notification Schedule: After Hours		Holiday Alarm Notification Schedules: Unassigned				
Input Type: Temperature Serial Number: Channel 1 Device Code: 3SS Input in use? Yes Expiration Date: 18-Aug-2017 Calibration Correction: +0.010		Pre-Alarm Standard Emergency				
		Alarm Min (C)		0.0	-5.00	0.0
		Alarm Max (C)		0.0	0.0	0.0
		Alarm Delay (minutes)		0	60.00	0
		Latching				
Sensor ID: fa00127e Sensor Type: CheckPoint		Alarm Limit In Use		x		
AA04873 - Lunch Room (Frz) (AA04873)	Reach-In	Lunch Room	global, Client Security, Device, and	Active		

System Configuration Report Example